

ASEPTIQUIK® G STC SERIES ASSEMBLY PROCEDURE

FLIP-CLICK-PULL Assembly Procedure

1

Steam flows from the process equipment through the Steam-Thru II in a steam trap creating a "steam on" sterile connection.

2

Unsnap and flip down the protective pull tab covers on each genderless AseptiQuik G connector half.

3

Align genderless AseptiQuik G couplings, push together until "CPC Click" confirmation.

4

Pull the snapped-together membrane pull tabs from the connector. Remove the Steam-Thru II lock sleeve.

5

Once the genderless AseptiQuik G assembly is complete and the steam trap has been closed, simply press the thumb latch to allow the valve to be pushed into the flow position. The "CPC Click" confirms the transition is complete.

6

Once the valve is locked into the flow position, you are ready for sterile fluid transfer to or from the process equipment.

7

After sterile media transfer is complete, simply press the thumb latch and pull the valve back into the steam position. The "CPC Click" confirms the transition is complete.

8

With the valve locked securely into the steam position, complete a second SIP cycle to "steam off" the connection.